

Please note: The concentration of the included pKanR has been changed from 1 ng/ μ l to 10pg/ μ l. Do not dilute the plasmid before performing the transformation positive control. Please contact Lucigen if you have any questions.

BAC-Optimized ReplicatorTM v2.0 Electrocompetent Cells

**IMPORTANT!
-80°C Storage Required
Immediately Upon Receipt**

Lucigen[®] Corporation
Advanced Products for Molecular Biology

2120 W. Greenview Drive
Middleton, WI 53562
Toll Free (888) 575-9695
Phone (608) 831-9011
FAX (608) 831-9012
E-mail lucigen@lucigen.com

www.lucigen.com

BAC-Optimized ReplicatorTM v2.0 Electrocompetent Cells

Contents

Components & Storage Conditions.....	3
BAC-Optimized Replicator v2.0 Electrocompetent Cells.....	3
Preparation for Transformation.....	4
Transformation/DNA Isolation.....	5
Media Recipes.....	6
Related Products.....	6

Notice of Limited Label License, CopyRight BAC, Patents, Warranties, Disclaimers and Trademarks

* The use of the CopyRight BAC Cloning Kits is covered under U.S. Patent No. 5,874,259 and other patents pending. This product is sold under license from the University of Wisconsin-Madison. Its use is limited solely to research purposes.

Copyright © 2001-2010 by Lucigen Corp. All rights reserved. UltraClone, CloneDirect, pEZ, pEZSeq, cSMART, and Replicator are trademarks of Lucigen Corp. Lucigen, BigEasy, CloneSmart, ClonePlex, CopyRight, DNATerminator, *E. coli*, PCRTerminator, and pSMART are registered trademarks of Lucigen Corp. HydroShear[®] and GeneMachines[®] are registered trademarks of GeneMachines Corporation (San Carlos, CA).

Lucigen's products are sold for research use only and are not to be used in humans or for medical diagnostics. Lucigen's liability with respect to any CloneSmart product is limited to the replacement of the product. No other warranties of any kind, expressed or implied, including without limitation, any implied fitness for any particular use, are provided by Lucigen. Lucigen is not liable for any direct, indirect, incidental or consequential damages arising out of or in connection with the use or inability to use any of its CloneSmart products.

Limited Label License

This product is the subject of U.S. Patent #5,874,259 and pending patent applications licensed to Lucigen Corporation. The consideration paid for this product grants a Limited License to use the product pursuant to the terms set forth in this Limited Label License. Academic, Not-for-Profit and For-Profit institutions acquire certain limited nontransferable rights with the purchase of this product (see below).

The purchase price of this product includes limited, nontransferable rights to use only the purchased amount of the product to perform CopyRight Cloning Technology and only as described in CopyRight and Replicator v2.0 Kit Instruction Manuals. **This limited license specifically excludes manufacture of the Replicator v2.0 competent cells.** Lucigen Corporation reserves all other rights; in particular, the purchaser of this product may not transfer or otherwise sell this product or its components or derivatives to a third party, and no rights are conveyed to the purchaser to use the product or its components or derivatives for commercial purposes. "Commercial purposes" includes any activity for which a party receives consideration and may include, but is not limited to, (1) use of the product or its components or derivatives in manufacturing, (2) use of the product or its components or derivatives for diagnostic purposes, (3) transfer or sale of vectors made with the product or components or derivatives of the product, (4) use of this product or materials made therefrom to provide a service, information, or data (e.g., DNA sequence) to a third party in return for a fee or other consideration, or (5) resale of the product or its components or derivatives, whether or not such product or its components or derivatives are resold for use in research. Academic, Not-for-Profit, and For-Profit institutions must obtain a separate license from Lucigen Corporation to use this product for any purpose other than those permitted above. It is the sole responsibility of the buyer to ensure that use of the product does not infringe that patent rights of third parties.

If the purchaser is not willing to accept these use limitations, Lucigen Corporation is willing to accept return of the product for a full refund. For information on obtaining a license, contact Lucigen Corporation, 2120 W. Greenview Dr., Middleton, WI 53562. Email: Lucigen@lucigen.com. Phone: 608-831-9011. Fax 608-831-9012.

BAC-Optimized Replicator™ v2.0 Electrocompetent Cells

Components & Storage Conditions

Lucigen's BAC-Optimized Replicator v2.0 Electrocompetent Cells yield $\geq 2 \times 10^{10}$ cfu/ μ g supercoiled pKanR DNA. The cells are shipped on dry ice in one container, along with supercoiled control pKanR DNA (10 μ g/ μ l), Arabinose Induction Solution, and Recovery Medium. BAC-Optimized Replicator v2.0 Cells are packaged in 50- μ l aliquots, sufficient for two transformations per tube. Please refer to the table below for catalog numbers.

Lucigen's BAC-Optimized Replicator v2.0 Electrocompetent Cells require storage at -80° C.

BAC-Optimized Replicator v2.0 Electrocompetent Cells Store at -80° C

	Catalog #	Reactions
BAC-Optimized Replicator v2.0 Electrocompetent Cells (DUOS) Store at -80° C	60210-1	12 (6 x 50 μ l)
	60210-2	24 (12 x 50 μ l)
Positive Control Plasmid (pKanR 10 μ g/ μ l) Store at -20° C to -80° C	----	(1 x 20 μ l)
Arabinose Induction Solution Store at -20° C to -80° C	----	1 ml
Recovery Medium Store at -20° C to -80° C	----	12 (1 x 12 ml)
	----	24 (2 x 12 ml)
	80026-1	96 (8 x 12 ml)
YT Agar (powder)	----	

BAC-Optimized Replicator v2.0 Electroporation Competent Cells

Lucigen's BAC-Optimized Replicator v2.0 Electrocompetent Cells are an *E. coli* strain that contains an inducible *trfA* gene, which is required for induction of replication from the *oriV* origin in BAC and fosmid clones. Most lab strains of *E. coli* do not contain a *trfA* gene, and thus will not support copy number amplification of such vectors.

BAC-Optimized Replicator v2.0 Cells are designed for high efficiency transformation by electroporation, yielding $\geq 2 \times 10^{10}$ cfu/ μ g pKanR DNA. These cells are ideal for cloning and propagation of BAC, fosmid, or plasmid clones. In addition to the *trfA* gene, they contain the *endA1* mutation for high yield and high quality plasmid DNA. They also contain the *mcr* and *mrr* mutations, stable cloning of methylated genomic DNA isolated directly from mammalian or plant cells. They do not contain the F plasmid. The *rpsL* mutation confers resistance to streptomycin.

BAC-Optimized Replicator v2.0 Genotype: F^{-} *mcrA* Δ (*mrr-hsdRMS-mcrBC*) *endA1* *recA1* ϕ 80d*lacZ* Δ M15 Δ *lacX74* *araD139* Δ (*ara,leu*)7697 *galJ galK rpsL nupG* (*attL* *araC-P*_{BAD}-*trfA250 bla attR*) λ^{-}

Note that Replicator v2.0 cells are resistant to ampicillin (*bla* gene).

The plasmid pKanR (kanamycin resistant) is provided as a control for transformation of Lucigen's Replicator v2.0 cells. Kanamycin is required for selection of pKanR transformants.

BAC-Optimized Replicator™ v2.0 Electrocompetent Cells

Preparation for Transformation

Large-insert DNA cloning and BAC library construction demand the highest transformation efficiency and recovery of the largest clones possible. Lucigen's BAC-Optimized Replicator v2.0 Electrocompetent Cells were developed exclusively for this purpose. These cells are provided in 100- μ l aliquots, sufficient for five transformation reactions. We also recommend use of the Cell Porator and Voltage Booster System from Whatman Biometra. Optimal settings for electroporation are listed below:

Table 2. Electroporation conditions for *E. cloni*® 10G BAC-Optimized Electrocompetent Cells

Optimal Setting (Recommended)	Alternate Conditions (May give lower efficiencies)
Cuvette gap: 1.5 mm Voltage: 358 V Capacitance: 330 μ F Impedance: Low ohms Charge rate: Fast Voltage Booster Resistance: 4000 ohms	Cuvette gap: 1.0 mm Voltage: 1800 V Capacitance: 25 μ F Impedance: 200 ohms
Whatman Biometra: Cell Porator and Voltage Booster System, Cat. # 11609013	Eppendorf: Model 2510; Bio-Rad: Gene Pulser II #165-2105; BTX: ECM630
1.5 mm Cuvettes: Whatman Biometra Cat.#11608031	1 mm Cuvettes: Eppendorf Model 4307-000-569, BTX Model 610, or BioRad Cat. #165-2089. Users have reported difficulties using <i>E. cloni</i> cells with Invitrogen cuvettes Cat. # 65-0030.

Optional transformation control reactions include electroporation with 1 μ l (10 pg) of supercoiled pKanR DNA.

To ensure successful transformation results, the following precautions must be taken:

- For best results, the BAC ligation reaction must NOT be purified or heat treated.
- The DNA sample to be used for electroporation must be dissolved in water or a buffer with low ionic strength, such as TE. The presence of salt in the electroporation sample leads to arcing at high voltage, resulting in the loss of the cells and DNA.
- Microcentrifuge tubes and electroporation cuvettes must be thoroughly pre-chilled on ice before use. Optimal results are obtained with the electroporator and cuvettes from Whatman Biometra (Cat.# 11609013 and #11608031, respectively). Successful transformation also can be achieved with alternate systems, although efficiency may be lower (see Table 2).
- The cells must be completely thawed **on ice** before use.
- For highest transformation efficiency, use the provided Recovery Medium to resuspend the cells after electroporation. Use of TB or SOC will result in lower transformation efficiencies.
- Use YT agar plus appropriate antibiotic for plating cells. Colony size will be small or variable on LB agar plates. YT Agar is provided to maximize colony size. Cells may be plated on LB or other common media—colonies will be noticeably smaller upon comparison but adequate for the vast majority of intents and purposes.

BAC-Optimized ReplicatorTM v2.0 Electrocompetent Cells

Transformation Protocol

1. Aliquot 1 ml of Recovery Medium into 17 mm x 100 mm sterile culture tubes at room temperature (one tube for each transformation reaction). Transformation efficiency may decrease with the use of other media.
2. Place electroporation cuvettes and microcentrifuge tubes on ice (one cuvette and one tube for each transformation reaction).
3. Remove *E. coli* cells from the -80°C freezer and place on wet ice until they thaw **completely** (10-20 minutes).
4. When cells are thawed, mix them by tapping gently. Add 20 µl of *E. coli* cells to the chilled microcentrifuge tube on ice.
5. If using a Lucigen BAC Cloning Kit, add 1 µl of the BAC ligation reaction directly to the 20 µl of cells on ice. Do NOT heat inactivate the ligation reaction. (Heat-inactivating the ligation reaction will reduce the quality of BAC cloning.) Stir briefly with pipet tip; **do not** pipet up and down to mix, which can introduce air bubbles and warm the cells. Using more than 1 µl of ligation mix may cause electrical arcing during electroporation.

For ligation reactions using other commercial kits, please refer to the manufacturer's instructions.
6. Carefully pipet the cell/DNA mixture into a chilled electroporation cuvette without introducing bubbles. Electroporate according to the conditions recommended on p.4.
7. After electroporation, quickly transfer the cells into the Recovery Medium in the culture tube at room temperature.
8. Place the tube in a shaking incubator at 250 rpm for 1 hour at 37°C.
9. Spread up to 100 µl of transformed cells on YT agar plates containing the appropriate antibiotic.
10. Incubate the plates overnight at 37°C.
11. Transformed clones can be further grown in TB or in any other rich culture medium.

DNA Isolation

Transformants are grown in TB medium. Stable inserts of 10-40 kb can be grown overnight with shaking at 37°C in the presence of 1X Arabinose Induction Solution. DNA minipreps can be performed by standard methods.

For BACs and unstable smaller inserts, it may be necessary to grow the cultures without induction to an OD₆₀₀ of 0.2-0.3. To reach this OD, it is convenient to grow the cultures overnight at 37°C without shaking. The following morning, dilute the cultures 2-10 fold, and grow at 37°C with shaking at 225 rpm for 30 minutes. For each ml of culture, add 1 µl of 1000 X Arabinose Induction Solution. Continue growth for 2-3 hours at 37°C with shaking at 225 rpm.

Prepare DNA minipreps according to standard protocols.

BAC-Optimized Replicator v2.0 Electrocompetent cells are *recA endA* deficient and will provide high quality plasmid DNA.

BAC-Optimized Replicator™ v2.0 Electrocompetent Cells

Media Recipes

TB Culture Medium for Growth of Transformants

Per liter:

- 11.8 g Bacto-tryptone
- 23.6 g yeast extract
- 9.4 g dipotassium hydrogen phosphate (anhydrous)
- 2.2 g potassium dihydrogen phosphate (anhydrous)
- 0.4 % glycerol

Add all components except glycerol to deionized water. Autoclave and cool to 55°C. Immediately before use, add 8 ml of filter-sterilized 50% glycerol plus appropriate antibiotics.

YT Agar Plates*

Add the YT Agar powder provided with the kit to 500 ml of deionized water. Autoclave and cool to 55°C. Add the appropriate filter-sterilized antibiotic to the cooled medium (e.g., 15 mg kanamycin for kanamycin-resistant transformants; 50 mg ampicillin or carbenicillin for ampicillin-resistant transformants). For blue/white screening, add 1.5 ml 100mM IPTG and 5 ml 2% X-gal to the molten agar at 55°C before pouring. Pour approximately 25 ml per petri plate.

Temperatures of >55°C may destroy the antibiotics. Do NOT add antibiotics to hot media!

Per liter YT is:

- 5 g yeast extract
- 8 g tryptone
- 5 g NaCl
- 15 g agar

Add deionized water to 1 liter. Adjust pH to 7.0 with NaOH.

*YT Agar is provided to maximize colony size. Cells may be plated on LB or other common media—colonies will be noticeably smaller upon comparison but adequate for the vast majority of intents and purposes.

Related Lucigen Products

- CopyRight® pSMART® VC Cloning Kits
- CopyRight pEZ BAC Cloning Kits
- BigEasy™ v2.0 Linear Cloning Kits
- CloneSmart® Blunt Cloning Kits
- DNATerminator® End Repair Kit
- UltraClone™ DNA Ligation & Transformation Kits
- CloneDirect™ Rapid Ligation Kit
- GC Cloning & Amplification Kits
- pEZSeq™ Blunt Cloning Kit
- cSMART™ cDNA Cloning Kit
- *E. coli*® Electrocompetent Cells